


Quick Start Card

To Crimp

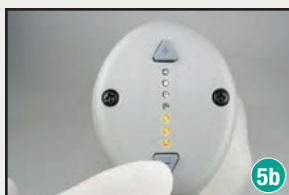

- 1 Attach a crimping jaw by screwing it on to the bottom of the unit. The crimping jaw can be identified by the downward facing arrow at the top of the jaw.
- 2 Apply power to the unit by plugging in the power supply cord at the top of the unit.
- 3 The compression of the Crimpenstein jaws will need to be adjusted based on the vials and seals that are being used. It is recommended that you start at the lowest crimping setting and increase the crimping compression as needed until the perfect crimp is achieved.
- 4 Rest the unit on top of the aluminum seal and press the trigger button to crimp.
- 5 Check the crimp and adjust the crimping jaw compression if necessary by pressing the "+" or "-" buttons on top of the unit.
 - a. To create a tighter crimp, press the "+" button on top of the unit. The lights will advance toward the red side of the display indicating a tighter crimp.
 - b. To create a looser crimp, press the "-" button on top of the unit. The lights will advance toward the yellow side of the display indicating a looser crimp.

If the Unit Jams Making a Crimp

Cause:

The cap has likely been over crimped and deformed.

Remedy:


1. Reset the unit by unplugging the power from the unit and then plugging the power back in. This will cause the unit to reverse and release the jaws.
2. If the above remedy fails to release the cap, gently press (do not bang) vial on a flat surface to force the jaws to release. Use the '-' button at the top of the unit to advance towards the yellow side of the display for a lighter crimp.
3. Use the '-' button at the top of the unit to adjust for a proper crimp.

Quick Start Card

To Decap

- 1 Attach a decapping jaw by screwing it on to the bottom of the unit. The decapping jaw can be identified by the upward facing arrow at the top of the jaw.
- 2 Apply power to the unit by plugging in the power supply cord at the top of the unit.
- 3 All the lights will illuminate when the unit is in decapping mode. Decapping requires the maximum power setting, and no adjustments are able to be made.
- 4 Rest the unit on top of the aluminum seal and press the trigger button to decap.

For more detailed information on crimping and decapping, please refer to the Crimpenstein Electric Crimper/Decapper instruction manual.

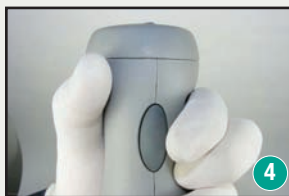

 **WHEATON** Science Products
a division of **WHEATON** Industries Inc.

1501 North 10th Street, Millville, NJ 08332-2038
800.225.1437 (U.S. & Canada) • 856.825.1100 • 856.825.1368 (F)
www.wheatonsci.com